




### TOPIC: Layout Save

The layout save feature allows the user to conveniently save multiple image files into one file for easier viewing and comparison.




- 1) Open up the PFV and go to the "FILE" tab
- 2) Select "Open File" and open the desired amount of files you wish to layout save. Ensure any post processing desired, such as: bit shift, gamma adjustments, HDR, etc., is completed as the layout save

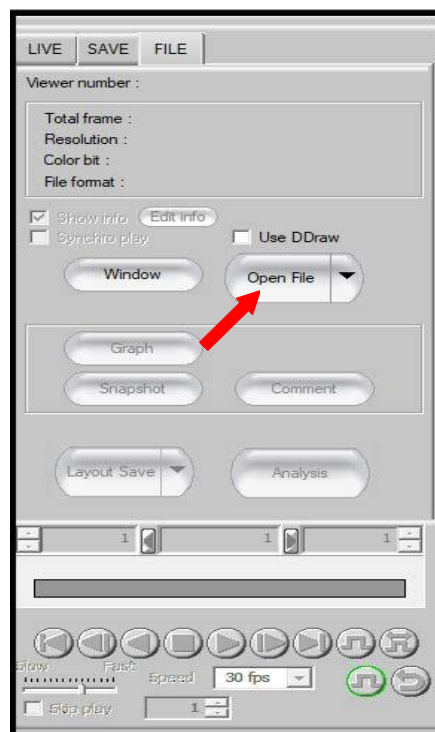


will only be 8-bit

3) Click on the drop down menu arrow and select "LayoutSave"

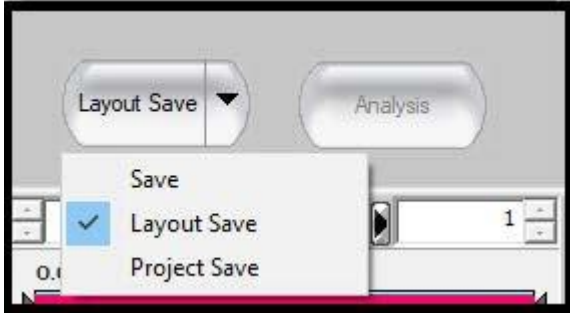


4) On the layout save menu, the user can select how they would like the image layout (In this example I selected 4 images and decided to use a 2x2 layout, but could also do 1x4 or 4x1). The order of images can also be manipulated by simply dragging the file name higher or lower in the "Target window(s)" section. The other options available are adjusting resolutions and saving desired information for each image.

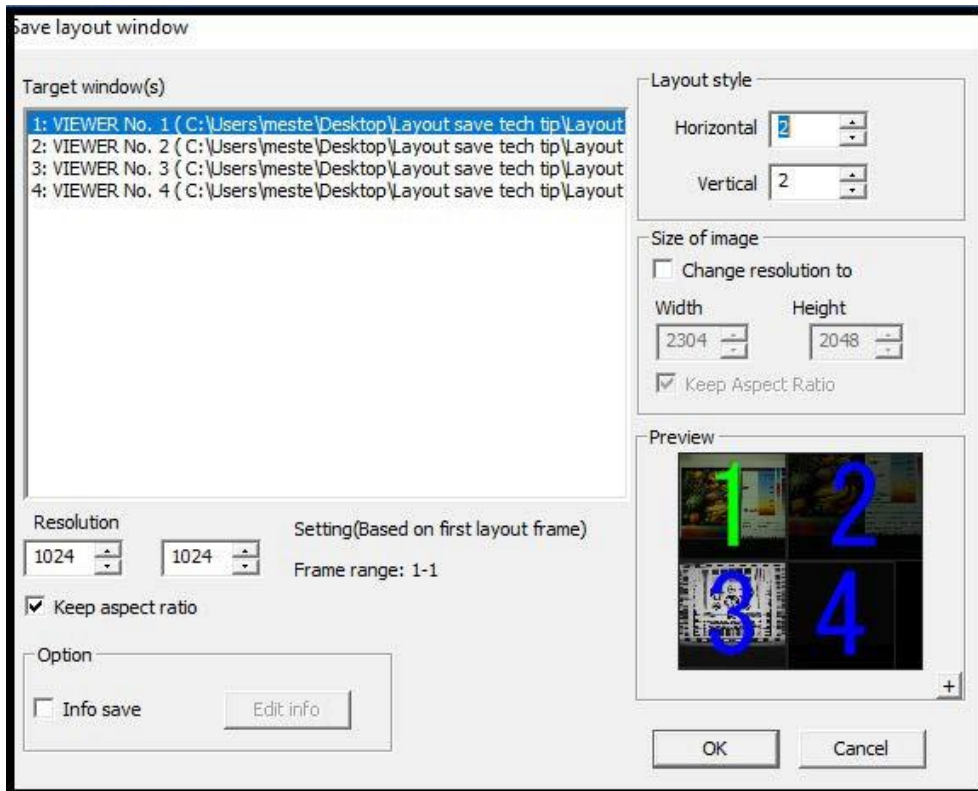

5) Select OK and choose a file name and file type (TIFF, BMP, JPEG, PNG, AVI, WMV) for your image. Please remember only 8 bits can be saved.

6) The layout save feature can also be used to save multiple videos. The videos must all contain the same number of frames and have the same trigger frame for layout save to work. A picture of the trigger frame adjustment button is pictured below. Select the button and move the red slider to your desired frame. Click on the down arrow to make that frame the trigger frame.

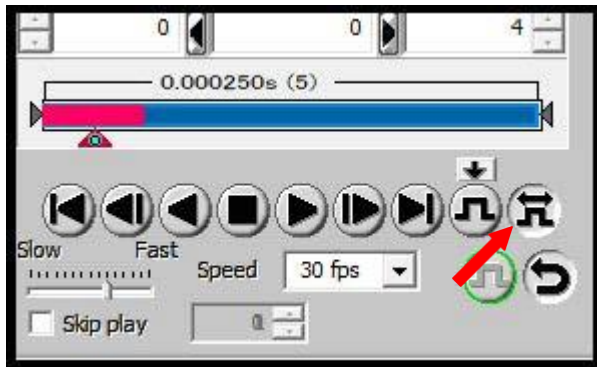

We hope this tech tip was helpful and if you have any questions or comments please contact us:  
Call: toll free at 800.585.2129 or 858.684.3555  
E-mail: [Mestela@photron.com](mailto:Mestela@photron.com) or [Skirksey@photron.com](mailto:Skirksey@photron.com)