



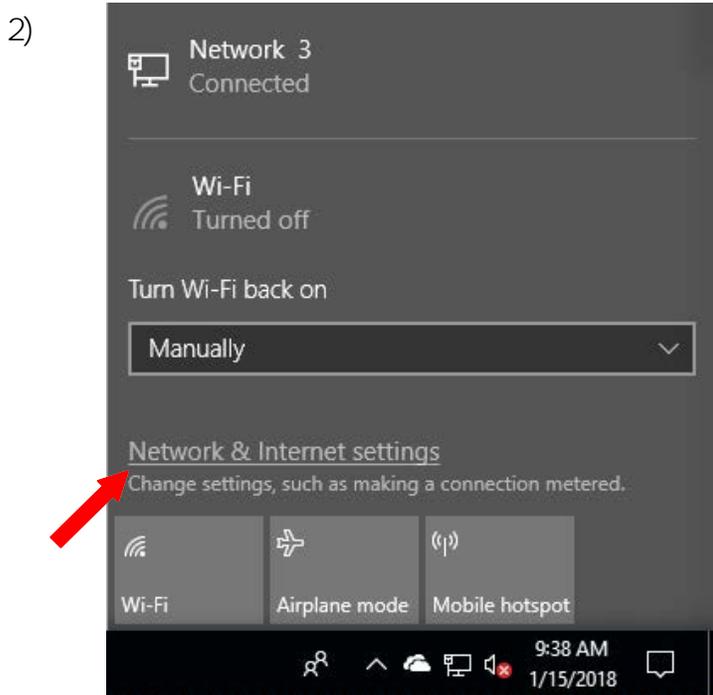

Timing mobile Intel network adapters for Windows 10

The Intel mobile i series of network adapters has issues communicating properly with Photron cameras if the firmware is out of date. Please ensure if you have an Intel adapter that starts with I you upgrade to the latest driver version.

To check your network adapter simply click on your ethernet or wifi icon in the bottom right of the task bar and select "Network and Internet Settings". The network settings window will open up and then the user can select "Change adapter options" to open up their list of adapters. Select the currently used adapter and right click and select properties. The name will be displayed in the box that says "Connect using:".

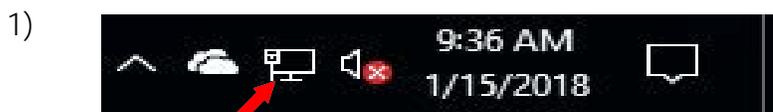



Status

3)

Network status



You're connected to the Internet

If you have a limited data plan, you can make this network a metered connection or change other properties.

[Change connection properties](#)

[Show available networks](#)

Change your network settings

-  **Change adapter options**
View network adapters and change connection settings.
-  **Sharing options**
For the networks you connect to, decide what you want to share.
-  **HomeGroup**
Set up a homegroup to share pictures, music, files, and printers with other PCs on your network.
-  **Network troubleshooter**
Diagnose and fix network problems.

